


# Human PBMC Isolation and Counting Using the Scepter™ 2.0 Handheld Automated Cell Counter

## Introduction

Peripheral blood mononuclear cells (PBMCs) are blood cells with round nuclei, such as monocytes and lymphocytes, with the lymphocyte population comprised of T cells, B cells, and natural killer (NK) cells. PBMCs are a critical component of the immune system, playing an integral role in the body's defense mechanisms. Cellular assays using PBMCs form the backbone of immune monitoring studies in clinical diagnostics and therapeutic design. Given that ineffective separation of lymphocytes from whole blood can significantly alter cellular responses and lead to unreliable results, it is essential to start every assay with a rapid, simple, and reliable method for PBMC isolation and subsequent quantitation.

Separation of PBMCs from whole blood is most commonly achieved via Ficoll® density gradient centrifugation.<sup>1-3</sup> Differential migration during centrifugation results in the separation of cell types into different layers (**Figure 1**). The bottom layer contains Ficoll-aggregated red blood cells. Immediately above this is a diffuse layer containing mostly granulocytes and unbound Ficoll®. Due to a slightly lower density, the lymphocytes (including the monocytic PBMC fraction) sediment at the interface between the Ficoll® and uppermost plasma/platelet layer. PBMCs are removed from the interface and subjected to multiple washes in PBS or cell medium to remove any residual Ficoll. Cell isolates are then ready for analysis or culture seeding.

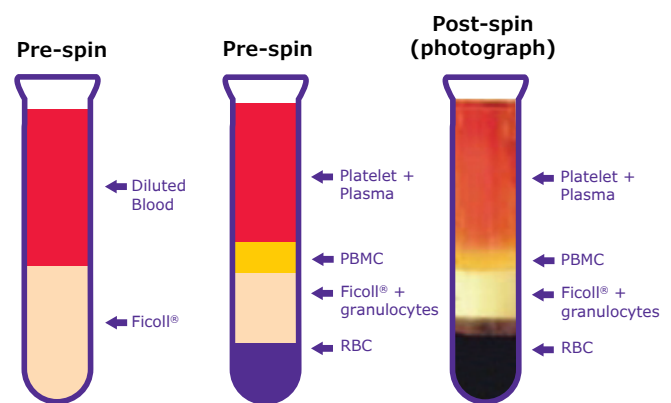

**Figure 1.** Diluted whole blood sample layering before and after Ficoll® density gradient centrifugation.

The Scepter™ cell counter combines the ease of an automated instrument and the accuracy of impedance-based particle detection using the Coulter principle in a handheld format. The instrument uses a combination of analog and digital hardware for sensing, signal processing, data storage, and graphical display. The precision-machined, consumable polymer sensor has a laser-drilled aperture in its cell sensing zone that enables the instrument to use the Coulter principle to discriminate cell diameter and volume at sub-micron and sub-picoliter resolution, respectively. Previous work has shown that, by using the 40 µm aperture sensor, the Scepter™ cell counter was able to accurately and precisely count a much broader range of cell types, including small cells (< 6 µm in diameter) such as PBMC and red blood cells (RBC).<sup>4</sup> This short report outlines a protocol for the isolation of PBMC from whole blood and subsequent sample analysis using the Scepter™ cell counter.

## Materials and Methods

This protocol is used to fractionate 10 mL of defibrinated or anticoagulant-treated peripheral blood\* (or buffy coat) using 15 mL Ficoll® separation medium in a 50 mL conical tube. A 15 mL tube can be used for fractionating smaller volumes of blood; however, for high PBMC yield with efficient RBC removal, it is important to maintain the same volumetric ratio of sample to Ficoll. Further, to ensure high viability of isolated cells, we recommend using only freshly isolated blood samples (< 24 hours post-collection).

\* Anticoagulants include: Heparin, EDTA, citrate, acid citrate dextrose (ACD), and citrate phosphate dextrose (CPD).

### Ficoll® density gradient separation of PBMC

1. Transfer 10 mL of blood from the collection vial to a 50 mL conical tube.
2. Add an equal volume of PBS (1X EmbryoMax® PBS, BSS-1006-A) and mix sample by repeated pipetting

**Note:** Diluting blood reduces the degree of RBC aggregation as well as freeing trapped PBMC. If left undiluted, trapped PBMC may sediment with erythrocytes, reducing yield.

3. Add 15 mL Ficoll® (GE Healthcare) to a second 50 mL tube.
4. Carefully layer the diluted blood over the Ficoll® by pipetting slowly and with minimal force.

**Note:** The diluted blood is added to the gradient by gently pipeting onto the separation medium with the tube held at an angle. To obtain good separation, it is paramount that clean separation of the blood and Ficoll® layers is maintained prior to centrifugation

5. Centrifuge without the brake applied at 400 g x 30 min at 18–24°C.

**Note:** Higher temperatures (37°C) enhance RBC aggregation and reduce yield, while lower temperatures (4°C) inhibit aggregation, decreasing purity. We recommend centrifuging at 18–24°C.

6. Carefully remove the tubes from the centrifuge so as not to disturb the layering.
7. Draw off the upper plasma layer being careful not to disturb the lower PBMC interface.
8. Remove the PBMC layer and transfer to a new 50 mL tube. The volume recovered should be approximately 10–12 mL.

**Note:** Attempt to collect the entire PMBC layer while avoiding the adjoining Ficoll(R) media and plasma layers. Recovery will result in contamination by granulocytes or platelets and plasma proteins, respectively.

9. Wash PBMC fraction using ~3 volumes of PBS  
Centrifuge at 100 g x 10 min at 18–24°C.
10. Decant the supernatant. Resuspend the pellet in 5 mL PBS. Fill PBS to 50 mL and repeat wash step.

11. Optional: The wash step can be repeated once more
12. Decant the supernatant and resuspend the cell pellet in appropriate volume of PBS (or media).

**Notes:** From healthy blood, PBMC yield ranges from 0.5 to 3 x 10<sup>6</sup> cells per mL blood. For 10 mL blood, resuspend pellet in 5 mL PBS for initial count. Purification of the PBMC population may be greatly enhanced by adding RosetteSep® Human Total Lymphocyte Enrichment Cocktail (StemCell Technologies Catalogue No. 15223) after this step.

13. Analyze samples using the Scepter™ cell counter and guava easyCyte™ flow cytometer.
14. Samples may require further dilution for accurate counting using the Scepter™ counter. The operating cell concentration range for the 40 µm aperture sensor = 5 x 10<sup>4</sup> – 1.5 x 10<sup>6</sup>.

### Scepter™ cell counting

The Scepter™ cell counter was used to count samples following the detailed on-screen instructions for each step of the counting process. Briefly, the user attaches a 40 µm sensor, depresses the plunger, submerges the sensor into the sample, then releases the plunger to draw 75 µL of cell suspension into the sensor. The Scepter™ cell counter detects each cell passing through the sensor's aperture, calculates cell concentration, and displays a size-based histogram as a function of cell diameter or volume on its screen. Scepter™ Software Pro was then used to upload files from the device and perform subsequent data analysis to determine the concentrations and relative cell frequencies for the lymphocyte and monocyte fractions.

### Guava easyCyte™ cell counting

Each 10 µL PBMC sample was diluted in 190 µL of PBS. Samples were then analyzed on a Guava easyCyte™ HT flow cytometer to determine the concentrations and relative cell frequencies for the lymphocyte and monocyte fractions.


### Cell surface staining and subset determination

For each sample, 100,000 PBMCs were resuspended in 100 µL PBS+0.1% BSA. To distinguish the discrete cell subsets present in PBMC, samples were stained with the following combination of fluorescently labeled antibodies: anti-CD3-PE (T-cells), anti-CD19-Alexa Fluor® 488 (B-cells), anti-CD16/CD56-APC (NK cells), and anti-CD14- PECy7 (monocytes) (eBioscience). Samples were incubated at room temperature for 20 minutes, washed with PBS, and resuspended in 200 µL PBS before acquisition. Samples were analyzed (3,000 cells/sample well) on a guava easyCyte™ HT flow cytometer using ExpressPro software.

## Results

Density gradient centrifugation using Ficoll® resulted in the characteristic separation of whole blood into 4 distinct layers: an uppermost plasma layer containing platelets, a thin PBMC band, a diffuse Ficoll® layer containing granulocytes, and the aggregated erythrocyte pellet. PBMC could be further subdivided into two main populations, the lymphocytes and monocytes.

Aliquots of each sample were stained with fluorescent antibody markers for characteristic PBMC phenotypes. The lymphocyte subset could be further subdivided into T cells (CD3+), B cells (CD19+), and NK cells (CD16/56+). Monocytes could be distinguished from total lymphocytes on the basis of differential CD14 expression. Samples were diluted and fractions analyzed using the flow cytometer and Scepter™ cell counter to determine total PBMC concentration.

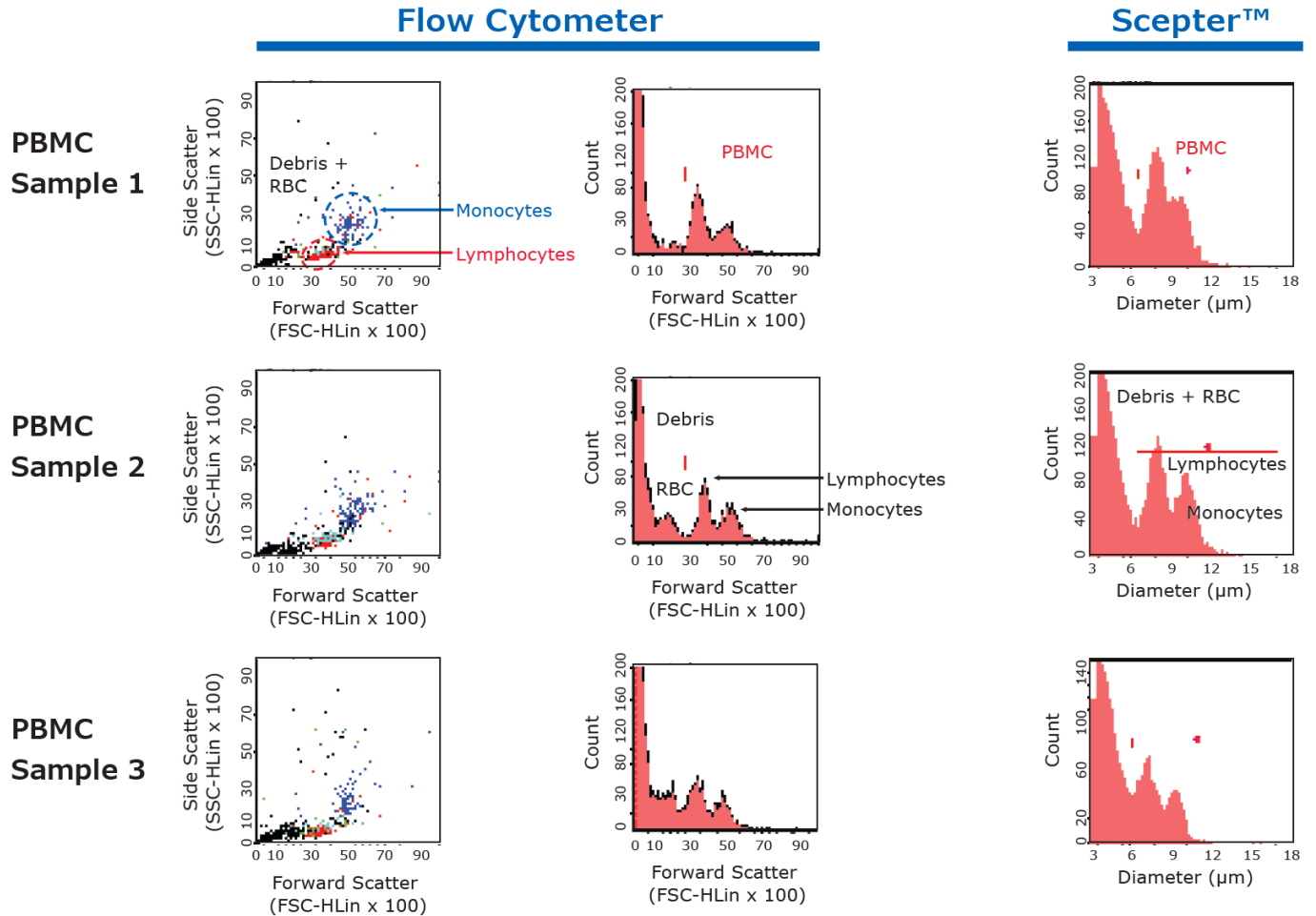

**Figure 2.**

Representative data comparing PBMC samples acquired on the flow cytometer and Scepter™ cell counter. Dot plots show debris

From the color dot plots in Figure 2, we discerned three event populations found within each sample: debris and red blood cells (Black), a lymphocyte subset (composed of T cells (Red), B cells (Aqua), NK cells (Green)), and a monocyte subset (Blue). While these subsets showed some overlap, each could be clearly defined by the size-based forward-scatter and diameter histogram plots of the flow cytometry and Scepter™ platforms, respectively. We observed somewhat greater peak distinction in the flow cytometry data than data from Scepter™ counting. Specifically, debris and contaminating erythrocytes each constituted a distinct subpopulation in flow cytometry derived histograms, but Scepter™ histograms showed only one peak corresponding to both fractions. Nine PBMC samples were analyzed (Table 1).

Across the nine samples, the average mean cell diameters were  $7.23 \pm 0.30 \mu\text{m}$  and  $10.02 \pm 0.20 \mu\text{m}$  for lymphocytes and monocytes, respectively. Resulting values are consistent with previously reported size ranges.<sup>5</sup> In addition, total PBMC concentrations were determined by three methods: Scepter™ diameter plot, flow cytometric forward scatter, and antibody staining (Figure 3). General agreement was observed amongst the different analytical techniques with values varying by <15% in all cases. Small differences in results may be due to subjectivity and user bias in the placement of gates defining the PBMC fraction.

### Relative frequency (percentage of total)

Test	Cell Fraction	Scepter™ cell countera	Forward Scatter <sup>b</sup>	Staining <sup>c</sup>
1	Lymphocyte	58	65	63
	Monocyte	42	35	37
2	Lymphocyte	68	72	71
	Monocyte	32	28	29
3	Lymphocyte	66	69	71
	Monocyte	34	31	29
4	Lymphocyte	62	67	64
	Monocyte	38	33	36
5	Lymphocyte	64	66	67
	Monocyte	36	34	33
6	Lymphocyte	62	58	60
	Monocyte	38	42	40
7	Lymphocyte	65	72	72
	Monocyte	35	28	28
8	Lymphocyte	59	61	61
	Monocyte	41	39	39
9	Lymphocyte	64	72	72
	Monocyte	36	28	28

**Table 1.** Lymphocyte and monocyte subset frequencies from nine individual PBMC samples. Aliquots from each sample were analyzed using the Guava easyCyte™ and Scepter™ instruments.

aValues were derived from the diameter histogram plot.


bValues were derived from the forward scatter histogram plot based on total events measured on guava easyCyte™ platform.

cStaining frequencies derived as follows:


%Lymphocytes = (%CD3+ T cells) + (%CD16/56+ NK cells) + (%CD19+ B cells)

%Monocytes = %CD14+ cells

PBMC Counts: Scepter™<sup>a</sup> cell counters vs. Flow Cytometry Histogram<sup>b</sup>

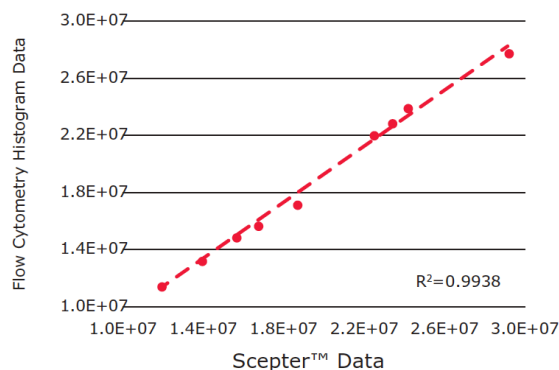

PBMC Counts: Scepter™<sup>a</sup> vs. Flow Cytometry Histogram<sup>b</sup>

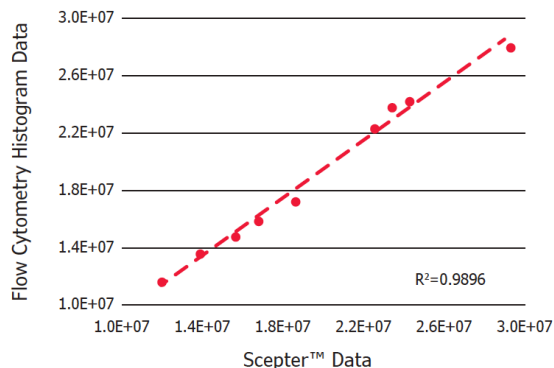

**Figure 3.** Correlation of PBMC concentrations measured using three different methods of analysis: flow cytometric cell counting, Scepter™ cell counting, and flow cytometry analysis of fluorescently labeled cells.

aValues were derived from the diameter histogram plot

bValues were derived from the forward scatter histogram plot based on total events

cPopulation counts derived as follows: %Lymphocytes = (%CD3+ T cells) + (%CD16/56+ NK cells) + (%CD19+ B cells); %Monocytes = %CD14+ cells

## Summary

PBMCs are widely used in both research and clinical laboratories. Separation of PBMCs from blood by centrifugation constitutes a critical first step for all downstream analyses. We have outlined the basic steps required for density-separation media based purification of mononuclear cells from whole blood and further shown that the Scepter™ cell counter can be used for accurate determination of resulting PBMC counts.

## Cited

1. Boyum, A. (1968) Isolation of mononuclear cells and granulocytes from human blood. Scand. J. Clin. Lab. Invest. 21:77-89.
2. Harris, R. and Ukaijiofo, E. (1970) Tissue typing using a routine one-step lymphocyte separation technique. British J. Haematol. 18:229- 35.
3. Bain, B. and Pshyk, K. (1972) Enhanced reactivity in mixed leukocyte cultures after separation of mononuclear cells on Ficoll-Hypaque. Transplantation Proceedings. 4:161-63.
4. Smith, J., and Ongena, K. The New Scepter™ 2.0 Cell Counter. Cellutions 2011 Vol. 1: p 19-22).
5. Daniels, V.G., Wheater, P.R., & Burkitt, H.G. (1979). Functional histology: A text and colour atlas. Edinburgh: Churchill Livingstone. ISBN 0-443-01657-7.

## General

- Prager, E., et al. (2001) Induction of Hyporesponsiveness and Impaired T Lymphocyte Activation by the CD31 Receptor: Ligand Pathway in T-Cells. J. Immunol. 166: 2364-2371.
- Smith, J., et al. The New Scepter™ 2.0 Cell Counter Enables the Analysis of a Wider Range of Cell Sizes and Types With High Precision. EMD Millipore Cellutions 2011 Vol. 1: p 19-22.

## To place an order or receive technical assistance

In Europe, please call Customer Service:

France: 0825 045 645	Spain: 901 516 645 Option 1
Germany: 069 86798021	Switzerland: 0848 645 645
Italy: 848 845 645	United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840

Or visit: [MerckMillipore.com/offices](https://www.MerckMillipore.com/offices)

For Technical Service visit: [MerckMillipore.com/techservice](https://www.MerckMillipore.com/techservice)

[MerckMillipore.com](https://www.MerckMillipore.com)

Merck KGaA  
Frankfurter Strasse 250  
64293 Darmstadt, Germany

