



## Product Information

### Trypsin-EDTA solution from porcine pancreas Cell Culture Tested

Product Number **T 4174**  
Storage Temperature -0 °C

#### Product Description

Enzyme Commission (EC) Number: 3.4.21.4  
Molecular Weight: 23.4 kDa<sup>1</sup>  
Extinction Coefficient:  $E^{1\%} = 15.0$  (280 nm)<sup>2</sup>  
pI: 10.2 - 10.8<sup>1,2</sup>  
Synonyms: Tryptase, Tryptar, Cocoonase,  
Parenzyme, Parenzymol

Trypsin Solution (5.0 g/l porcine trypsin and 2.0 g/l  
EDTA•4Na in 0.9% sodium chloride, 10X, cell culture  
tested)

Trypsin consists of a single chain polypeptide of  
223 amino acid residues. Trypsin is produced by the  
removal of the N-terminal hexapeptide from  
trypsinogen which is cleaved at the Lys<sup>6</sup> - Ile<sup>7</sup> peptide  
bond. The amino acid sequence of trypsin is cross-  
linked by 6 disulfide bridges. This native form of  
trypsin is referred to as  $\beta$ -trypsin. Autolysis of  $\beta$ -trypsin  
(which is cleaved at Lys<sup>131</sup> - Ser<sup>132</sup>) results in  $\alpha$ -trypsin  
which is held together by disulfide bridges. Trypsin is a  
member of the serine protease family. The active site  
amino acid residues of trypsin include His<sup>46</sup> and  
Ser<sup>183</sup>.<sup>1,3</sup>

Trypsin will cleave peptides on the C-terminal side of  
lysine and arginine amino acid residues. The rate of  
hydrolysis is slower if an acidic residue is on either  
side of the cleavage site and no cleavage occurs if a  
proline residue is on the carboxyl side of the cleavage  
site. The pH optimum of trypsin is 7 - 9.<sup>2</sup> Trypsin will  
also hydrolyze ester and amide linkages of synthetic  
derivatives of amino acids such as: benzoyl L-arginine  
ethyl ester (BAEE), p-toluenesulfonyl-L-arginine  
methyl ester (TAME), tosyl-L-arginine methyl ester,  
N $\alpha$ -benzoyl-L-arginine p-nitroanilide (BAPNA),  
L-lysyl-p-nitroanilide, and benzoyl-L-arginamide.<sup>2,3,4,5</sup>  
Assuming the pH and temperature are the same and  
using a molar extinction coefficient of 808 at 254 nm  
for BAEE, the following conversions are valid:

1 BAEE  $\mu$ M Unit = 200 BAEE Units  
1 TAME  $\mu$ M Unit = 0.27 BAEE  $\mu$ M Units  
1 BAEE  $\mu$ M Unit = 3.64 TAME Units  
1 TAME  $\mu$ M Unit = 55 BAEE A<sub>253</sub> Units  
1 BAEE A<sub>253</sub> Unit = 0.018 TAME  $\mu$ M Unit  
1 TAME  $\mu$ M Unit = 180 TAME A<sub>247</sub> Units  
1 TAME A<sub>247</sub> Unit = 0.33 BAEE Units  
A USP Unit =  $\Delta A_{253}$  of 0.003 per minute  
1 NF Unit = 3.3 A<sub>253</sub> BAEE Units.<sup>9</sup>

Note: These activity conversions were determined  
using bovine trypsin; however, they are thought to be  
similar for porcine trypsin.

Serine protease inhibitors that will inhibit trypsin  
include DFP (diisopropyl fluorophosphate),  
TLCK (N $\alpha$ -p-tosyl-L-lysine chloromethyl ketone),  
PMSF (phenylmethanesulfonyl fluoride),  
APMSF(4-amidinophenylmethanesulfonyl fluoride),  
AEBSEF (4-(2-aminoethyl)benzenesulfonyl fluoride),  
aprotinin, leupeptin,  $\alpha_2$ -macroglobulin,  $\alpha_1$ -antitrypsin,  
p-aminobenzamidine, benzamidine (reversible),  
soybean trypsin inhibitor, lima bean inhibitor, bovine  
pancreas trypsin inhibitor, chicken egg white inhibitor,  
and turkey egg white inhibitor.<sup>1,7</sup>

#### Precautions and Disclaimer

For Laboratory Use Only. Not for drug, household or  
other uses.

#### Storage/Stability

This product is stored frozen.

Trypsin retains most of its activity in 2.0 M urea,  
2.0 M guanidine HCl, or 0.1% (w/v) SDS.<sup>8</sup> Trypsin is  
reversibly denatured at high pH (above 11), by  
precipitation with TCA, or by high concentrations of  
urea (greater than 6.5 M).<sup>3</sup> In order to abolish all  
trypsin activity, heating at 100 °C in 1% (w/v) SDS for  
5 minutes is required.<sup>9</sup>

## Procedure

Trypsin may be used to remove adherent cells from a culture surface. Cells are most commonly removed from the culture substrate by treatment with trypsin, or trypsin•EDTA. Trypsin 1X solutions can range from 0.025% to 0.5%. The reasons for the range of concentrations are as follows: (1) Differences in trypsin activity or potency; (2) Different incubation times; (3) Different cell lines.

Incubating cells with too high a trypsin concentration for too long a time period will damage cell membranes and kill the cells. If unsure about the concentration of trypsin to use, use a low concentration. There can be lot-to-lot variation in dissociation times which is to be expected since the enzymatic activity of each lot will differ. If trypsin is being solubilized or diluted from a concentrated solution, it is important to use a buffered salt solution that contains no  $\text{Ca}^{2+}$  or  $\text{Mg}^{2+}$ , such as Hank's Balanced Salt Solution, Modified (Product No. H 9394). Adjust the pH of trypsin solution to 7.4-7.6.

- 1) Remove medium from culture vessel by aspiration and wash the monolayer with  $\text{Ca}^{+2}$  and  $\text{Mg}^{+2}$  - free salt solution to remove all traces of serum. Remove salt solution by aspiration.
- 2) Dispense enough trypsin or trypsin•EDTA solution into culture vessel(s) to completely cover the monolayer of cells and place in 37 °C incubator for approximately 2 minutes.
- 3) Remove the trypsin or trypsin•EDTA solution by aspiration and return closed culture vessel(s) to incubator. The coated cells are allowed to incubate until cells detach from the surface. Progress can be checked by examination with an inverted microscope.

Note: The time required to remove cells from the culture surface is dependent on cell type, population density, serum concentration in the growth medium, potency of trypsin and time since last subculture. Trypsin causes cellular damage and time of exposure should be kept to a minimum.

- 4) When trypsinization process is complete the cells will be in suspension and appear rounded.
- 5) It is advisable to add serum or medium containing serum to the cell suspension as soon as possible to inhibit further tryptic activity which may damage cells. Soybean trypsin inhibitor (Product No. T 9008) can also be added at an equimolar concentration to inhibit the trypsin that is present. Soybean trypsin inhibitor is used when culturing in serum-free conditions.
- 6) Cells can be resuspended by gently pipetting the cell suspension to break up the clumps. Further dilution can be made, if required, for cell counts and/or subculturing.

## References

1. Walsh, K. A., Trypsinogens and trypsins of various species. *Meth. Enzymol.*, **19**, 41-63 (1970)
2. Buck, F. F., et al., On the mechanism of enzyme action. LXXII. Studies on trypsins from beef, sheep, and pig pancreas. *Arch. Biochem. Biophys.*, **97**, 417-424 (1962).
3. Keil, B., in *The Enzymes*, 3rd ed., Vol. III, Boyer, P. D., Academic Press (New York, NY: 1971), pp. 250-275.
4. Burdon, R. H., et al., in *Laboratory Techniques in Biochemistry and Molecular Biology*, Vol. 9, 2nd ed., Allen, G., ed., Elsevier/North (New York, NY: 1989), pp. 73-104.
5. *Enzyme Handbook*, Vol. II, Barman, T. E., Springer-Verlag (New York, NY: 1969), pp. 618-619.
6. USP, Vol. 23, p. 1611.
7. *Proteolytic Enzymes, A Practical Approach*, Beynon, R. J., ed., IRL Press (New York, NY: 1989), p. 240.
8. *Methods of Molecular Biology*, Vol. 3, Smith, B. J., Humana Press, (New Jersey, 1988), pp. 57-69.
9. Porter, W. H., and Preston, J. L., Retention of trypsin and chymotrypsin proteolytic activity in sodium dodecyl sulfate solutions. *Anal. Biochem.*, **66**, 69-77 (1975).

TMG/RXR 8/03

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.